

$(\text{NH}_4)_2\text{SO}_4$

行业： 化学

进料尺寸：

最终精度： <400 μm

样品量： 85g

研磨建议：

Result 1



PLANETARY MONO MILL PULVERISETTE 6 CLASSIC LINE

600 rpm

500 ml ZrO₂
25x20 mm ZrO₂

Feed quantity: 250 g
Grinding time: 3 min

Result 2



MORTAR GRINDER PULVERISETTE 2

grinding set made of agate
pressure: 10 kg

Feed quantity: 85 g
Grinding time: 10 min
Final fineness: < app. 400 μm , sticks*1
Comments: *1: as the sample is hygroscopic it already sticks after a short grinding time on grinding bowl and grinding balls. A longer dry grinding is not possible.