

### Mycoplasma Off™

**Surface disinfectant for the complete elimination of mycoplasma from any cell culture work area and microbiology lab.**

#### Description

Mycoplasma Off™ is a simple solution for cleaning and disinfecting a variety of laboratory surfaces at the same time. There's no need for multiple products when you have the convenience of the easy-to-use Mycoplasma Off™ product line. Then using a simple sponge or paper towel, you may spread germs from one surface to another. Thus, germs are, or may be, propagated rather than wiped off.

Mycoplasma Off™ kills mycoplasma within seconds. It is also effective against common bacteria such as *Staphylococcus aureus*, *Streptococcus faecalis*, *Escherichia coli*, *Proteus mirabilis*, *Pseudomonas aeruginosa*, *Candida albicans*, *Salmonella enterica* and TBC. It kills a wide range of viruses including those causing colds and flu, HBV/HIV, Papova, Adeno, Polio and Vaccinia viridae, according to the requirements of CEN TC 216.

#### Protection and Precaution Information

Do not use for alcohol-sensitive surfaces. Pre-test the surface for color fastness in an inconspicuous area. Wear disposable gloves and safety glasses when using the product. Ensure adequate ventilation after usage. For research use only.



#### Instruction for Use

For disinfecting and cleaning of surfaces and laboratory equipment. Spray Mycoplasma Off™ directly onto the surface and dry with paper towels.

For decontamination of **electronic laboratory devices**, we recommend our Mycoplasma Off™ Wipes (Cat. 15-1001 and 15-5001).

**Composition:** 32.5 % 1-propanol, 18 % ethanol, 0.1 % glutaraldehyde, 0.05 % Mynox®

**Catalog numbers:** 15-1000 (1000 ml Spray).  
15-5000 (5 x 1000 ml Refill)

Germany:

**Minerva Biolabs GmbH**

Schkopauer Ring 13 | 12681 Berlin | Germany

Tel.: +49-30-2000437-0 | Fax: +49-30-2000437-9

Email: [order@minerva-biolabs.com](mailto:order@minerva-biolabs.com)

Internet: [www.minerva-biolabs.com](http://www.minerva-biolabs.com)

USA / Canada

**Minerva Biolabs Inc.**

1 Jill Ct., Building 16, Unit 10

Hillsborough | NJ 08844 | USA

Phone: 1-908-524-4661

Internet: [www.minervabiolabs.us](http://www.minervabiolabs.us)